



**Advantech Manufacturing, Inc**

2450 S. Commerce Drive  
New Berlin, WI 53151

Toll Free: 800.511.2096  
Tel: 262.786.1600  
Fax: 262.786.5074

[www.advantechmfg.com](http://www.advantechmfg.com)  
[sales@advantechmfg.com](mailto:sales@advantechmfg.com)

## **CARE AND CLEANING FOR METAL FRAMED WIRE CLOTH SIEVES**

*(For care and handling of electroformed nickel precision sieves,  
please refer to: CARE AND CLEANING OF PRECISION SIEVES.)*

- Clean sieves from **BOTTOM** side with a brush. Use extreme care to avoid damage.
- Never use sharp objects to clean your wire cloth sieves.
- Wash in a mild solution of detergent and warm water. For a more thorough cleaning, some ultrasonic\* cleaners may be used. *(\*Caution: Refer to your ultrasonic manufacturer instructions prior to use as some devices may cause screen damage.)*
- **CAUTION:** Do not expose sieves to temperatures above 261° Fahrenheit as solder will begin to soften.
- Do not use forced air to dry #100 or finer wire cloth mesh sieves. This may result in screen damage.
- Store equipment in a clean, dry environment when not in use.