



**Advantech Manufacturing, Inc**

2450 S. Commerce Drive  
New Berlin, WI 53151

Toll Free: 800.511.2096

Tel: 262.786.1600

Fax: 262.786.5074

[www.advantechmfg.com](http://www.advantechmfg.com)  
[sales@advantechmfg.com](mailto:sales@advantechmfg.com)



**CARE AND CLEANING FOR PRECISION SIEVES\***

**CAUTION – Electroformed Nickel Mesh  
Sieves are extremely fragile!!!**

Precision sieves incorporate an unsupported electroformed nickel mesh, which are very fragile and require special handling. Long sieve life can be obtained by observing the following guidelines:

- **Never** touch the sieve mesh with your fingers or objects other than your test media. Acids in perspiration will permanently stain or damage your screen.
- **Never** use a brush or forced air to clean precision mesh.
- **Never** expose precision sieves or mesh to radiant energy or temperatures above 120°F.
- Wash in a mild solution of detergent and warm water (70°F to 85°F). When cleaning, hold the sieve in a vertical position and gently rinse with distilled water only. Allow to air dry.
- Always store precision sieves in the original protective box, in a clean, dry environment.

***\* Sieve or mesh damaged caused by  
improper use, cleaning, or handling will  
void the manufacturer's warranty.***