



Advantech Manufacturing, Inc

2450 S. Commerce Drive
New Berlin, WI 53151

Toll Free: 800.511.2096

Tel: 262.786.1600

Fax: 262.786.5074

www.advantechmfg.com
sales@advantechmfg.com

CARE AND CLEANING FOR SONIC SIFTER WIRE CLOTH SIEVES

*(For care and handling of electroformed nickel precision sieves,
please refer to: CARE AND CLEANING OF PRECISION SIEVES.)*

- Clean sieves from **BOTTOM** side with a brush. Use extreme care to avoid damage.
- Never use sharp objects to clean your wire cloth sieves.
- Wash in a mild solution of detergent and warm water.
- Mating parts of sieves must be clean at all times to ensure a tight seal during sieving operations.
- **CAUTION:** Do not expose sieves to temperatures above 120° Fahrenheit.
- Do not use acids or cleaning agents that may damage the acrylic frame.
- Do not use forced air to dry #100 or finer wire cloth mesh sieves. This may result in screen damage.
- Store equipment in a clean, dry environment when not in use.